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Paraffin Sides Experiment Report

1. Experimental Equipment and Reagents

1.1 Experimental equipment

Name	Manufactor	Model
Dehydrator	DIAPATH	Donatello
Embedding machine	Wuhan Junjie Electronics Co., Ltd	JB-P5
Pathology slicer	Shanghai Leica Instrument Co., Ltd	RM2016
Frozen platform	Wuhan Junjie Electronics Co., Ltd	JB-L5
Tissue spreader	Zhejiang Kehua Instrument Co., Ltd	KD-P
Oven	Tianjin Laibo Rui Instrument Equipment Co., Ltd	GFL-230
Slides	Servicebio	

1.2 Main experimental reagents

Reagent	Manufactor	Article number
Xylene	Sinopharm Group Chemical Reagent Co. Ltd	10023418
Ethanol	Sinopharm Group Chemical Reagent Co. Ltd	100092683

2. Paraffin section making

2.1 The fresh tissue was fixed with fixed liquid for more than 24 hours. Remove the tissue from the fixed liquid and trim the target tissue with a scalpel in the ventilation cupboard, and put the trimmed tissue and the label in the dehydration box.

2.2 Dehydration and wax leaching: put the dehydration box into the dehydrator in order to dehydrate with gradient alcohol. 75% alcohol 4 hours, 85% alcohol 2 hours, 90% alcohol 2 hours, 95% alcohol 1 hour, anhydrous ethanol I 30 min, anhydrous ethanol II 30 min, alcohol benzene 5~10 min, xylene II 5~10 min, 65°C melting paraffin I 1h, 65 °C melting paraffin II 1h, 65°C melting paraffin III 1 hour.

2.3 The wax-soaked tissue is embedded in the embedding machine. First, put the melted wax into the embedding frame, and before the wax solidifies, remove the tissue from the dewatering box and put it into the embedding frame according to the requirements of the embedding surface and affix the corresponding label. Cool at -20 °C freezing table, and after the wax is solidified, the wax block is removed from the embedded frame and repaired.

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Wuhan Servicebio Technology Co., Ltd.

2.4 Place the trimmed wax block cool at 20° freezing table, slice the modified tissue chip wax block on the paraffin slicer, the slice thickness is 4μm. The tissue is flattened when the slice floats on the 40 °C warm water of the spreading machine, and the tissue is picked up by the glass slides and baked in the oven at 60 °C. After the water-baked dried wax is melted, it is taken out and stored at room temperature.