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Add: 5th Floor, C22 Building, Biopark, No. 388 Gaoxin 2nd Road, East Lake High-tech Developing Zone,
Wuhan, Hubei, China, 430079 Website: <http://www.servicebio.com> Tel: 027-5111-3188

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Acid-fast Staining Experimental Report

1. Lab Equipment and Reagents

1.1 Lab equipment

Items	Manufacturer	Model
Dehydrator	DIAPATH	Donatello
embedding machine	Wuhan Junjie Electronics Co., Ltd.	JB-P5
Pathology microtome	Shanghai Leica Instruments Co., Ltd.	RM2016
Frozen platform	Wuhan Junjie Electronics Co., Ltd.	JB-L5
Water Bath-Slide Drier	Zhejiang Jinhua Kedi Instrumental Equipment CO.,LTD	KD-P
Laboratory oven	Tianjin Labotery Instrument Equipment Co., Ltd.	GFL-230
Microscope slide	Servicebio	G6004
Upright optical microscope	Nikon Japan	Nikon Eclipse E100
Imaging system	Nikon Japan	NIKON DS-U3

1.2 Chemical reagents

Items	Manufacturer	Model
Absolute alcohol	Sinopharm Chemical Reagent Co., Ltd.	100092683
Xylene	Sinopharm Chemical Reagent Co., Ltd.	10023418
Acid-fast staining kit	Servicebio	G1047
Differentiation solution	.Servicebio	G1039
	Servicebio	G1040
Neutral balsam	Sinopharm Chemical Reagent Co., Ltd.	10004160

2. Experimental Steps

Preparation of Acid-fast staining solution: Acid-fast staining solution A 10ml + Acid-fast staining solution B 90ml, well mixed.

Preparation of acid-fast differentiation solution: take out 10ml of concentrated hydrochloric acid and dilute to 1000ml with absolute ethanol.

- 2.1 Paraffin section deparaffinize and rehydrate: put the slides into xylene I 20minutes-xylene II 20 minutes-absolute ethanol I 5 min-absolute ethanol II 5 min-75% alcohol for 5 min, then tap water washing.
- 2.2 Acid-fast staining: put the slides into the Acid-fast staining solution stain for 30 minutes and running water washing;
- 2.3 Background differentiation: acid-fast differentiation solution rapidly differentiates for 2s. Terminate differentiation by water washing, which can be repeated. By the time of microscope examination, the acid-fast bacteria are purple-red, and the background is basically colorless;
- 2.4 Sulphuric acid differentiation: If the background is deep and the hydrochloric acid differentiation does not fall, put the slides into acid-red staining C for rapid differentiation for 1s, tap water washing (repeatedly rinse until no sulfuric acid remains on the slide), and repeat operations until the positive bacteria are purple-red , the background is colorless.
- 2.5 Acid-fast solution D staining: put the slides into acid-fast solution D stain for 30s, tap water washing, differentiate for 2s, tap water washing, return blue solution to blue for 3-5s, rinse with running water.
- 2.6 Dehydration and sealing: put the slides into absolute ethanol I 5min-anhydrous ethanol II 5min-anhydrous ethanol III 5min-xylene I 5min-xylene II 5min -transparent, neutral balsam sealing.
- 2.7 Microscope examination, images collection and analysis.

3. Results

The acid-fast bacteria are purple-red filaments, and the nucleus is light blue.

4. Note

- 4.1 Pay attention to the degree of differentiation, over-differentiate, and differentiate the purple color of acid-fast bacteria., insufficient differentiation, the background is red.
- 4.2 Light staining during hematoxylin counterstaining process, should not affect the observation of acid-fast bacteria.