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Add: 5th Floor, C22 Building, Biopark, No. 388 Gaoxin 2nd Road, East Lake High-tech Developing Zone, Wuhan, Hubei, China, 430079 Website: <http://www.servicebio.com> Tel: 027-5111-3188

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## Copper Red Acid Staining Experimental Report

### 1. Lab Equipment and Reagents

#### 1.1 Lab equipment

| Items                      | Manufacturer                                           | Model              |
|----------------------------|--------------------------------------------------------|--------------------|
| Dehydrator                 | DIAPATH                                                | Donatello          |
| embedding machine          | Wuhan Junjie Electronics Co., Ltd.                     | JB-P5              |
| Pathology microtome        | Shanghai Leica Instruments Co., Ltd.                   | RM2016             |
| Frozen platform            | Wuhan Junjie Electronics Co., Ltd.                     | JB-L5              |
| Water Bath-Slide Drier     | Zhejiang Jinhua Kedi Instrumental<br>Equipment CO.,LTD | KD-P               |
| Laboratory oven            | Tianjin Labotery Instrument Equipment<br>Co., Ltd.     | GFL-230            |
| Microscope slide           | Servicebio                                             |                    |
| Upright optical microscope | Nikon Japan                                            | Nikon Eclipse E100 |
| Imaging system             | Nikon Japan                                            | NIKON DS-U3        |

#### 1.2 Chemical reagents

| Items            | Manufacturer                         | Model     |
|------------------|--------------------------------------|-----------|
| Absolute alcohol | Sinopharm Chemical Reagent Co., Ltd. | 100092683 |
| Xylene           | Sinopharm Chemical Reagent Co., Ltd. | 10023418  |
| Dithio oxamide   | Sigma                                | 379387    |
| Acetic acid      | Sinopharm Chemical Reagent Co., Ltd. | 10018818  |
| Nuclear fast red | Servicebio                           | G1035     |
| Neutral balsam   | Sinopharm Chemical Reagent Co., Ltd. | 10004160  |

### 2. Reagent Preparation

0.1% Dithio oxamide ethanol solution: Dithio oxamide 10mg + absolute ethanol 10ml

10% acetic acid: acetic acid 10g + ultra-pure water 100ml

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Add: 5th Floor, C22 Building, Biopark, No. 388 Gaoxin 2nd Road, East Lake High-tech Developing Zone,  
Wuhan, Hubei, China, 430079 Website: <http://www.servicebio.com> Tel: 027-5111-3188

3.1 Paraffin section deparaffinization and rehydration: put the slides into xylene I 20minutes-xylene II 20 minutes-absolute ethanol I 5 min-absolute ethanol II 5 min-75% alcohol for 5 min, then tap water washing and distilled water washing for 3 times.

3.2 Copper red acid staining: 0.1% Dithio oxamide ethanol solution: 10% sodium acetate mix at a ratio of 1:20, and put the slides into the mixed solution and covered in a 37°C water bath for about 3 days. Put the slides into 70% ethanol twice for 10 minutes each time, and absolute ethanol washing for twice, 3 hours each time;

3.3 Nuclear staining: distilled water washing for 3 times, wash out ethanol, nuclear fast red counterstaining for 3min, then tap water washing;

3.4 Dehydration and sealing: dehydrate with absolute ethanol in 3 cylinders, transparent with n-butanol for 5min, and xylene for 5min, neutral balsam sealing.

3.5 Microscope examination, image collection and analysis.

#### 4. Results

The copper salt deposition are dark green and black, and the nucleus is red.

#### 5. Notes

5.1 0.1% Dithio oxamide ethanol solution should be stored at room temperature( lightproof);

5.2 The copper red acid mixed solution should be ready for use and discard after usage.