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TTC Staining Experiment Report

I. Experimental Equipment and Reagents

1. Experimental equipment

Equipment name	Manufacturer	Model No.
Thermostatic water bath	Jiangyan Wuli Medical Instrument Factory Co., Ltd.	TL-42OD
Specimen bottle	Servicebio	
Leica blade	Shanghai Leica Instrument Co., Ltd.	LEICA 819
Dissecting forceps		
Camera	Canon	70D

2. Main experimental reagents

Reagent name	Manufacturer	Item No.
TTC Staining solution	Servicebio	G1017
Fixative	Servicebio	G1101

II. Preparation and staining steps of fresh tissue sections

1. Draw materials: After the animal is anesthetized, immediately take fresh tissues (such as brain and heart), then put the tissues without fixation directly into OCT embedding medium and quick-freeze at -20 °C for 30min. If not used urgently, the tissues can be temporarily cryopreserved at -80 °C.

2. Tissue block sectioning: Thaw the frozen tissue slightly and manually slice with a new Leica blade. The thickness of 2-3mm is suitable, and the cut surface should be neat.

3. TTC staining: Put the sections into the 37°C preheated TTC incubation solution, incubate in a 37°C water bath in the dark for 30 min, shake the tissues gently at 10 min intervals to make it evenly contact with the staining solution, and Observe the staining results.

4. Photographing: After the tissues are in color, add into fixative to stop the reaction, then preserve the tissue in dark, and take photo with camera.

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Add: 5th Floor, C22 Building, Biopark, No. 388 Gaoxin 2nd Road, East Lake High-tech Developing Zone, Wuhan, Hubei, China, 430079 Website: <http://www.servicebio.com> Tel: 027-5111-3188

The normal tissue area is red, and the infarct area is pale.

IV. Precautions:

1. The fresher the specimen, the better, do not fix, the blade should be thin and sharp, and the time of quick freezing in refrigerator should be suitable for sectioning smoothly, keep the sections flat.
2. The stained specimens should be stored in fixative and protected from light, then the staining results can be kept for several months, they are easily faded in the sun.
3. The TTC staining solution can be stored for one month in the dark at 4°C, return to normal temperature before use.

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Wuhan Servicebio Technology Co., Ltd.