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Experimental Report on Decalcification and Paraffin Embedding of Bone Tissue

1. Experimental equipments and reagents

1.1 Experimental equipments

Name	Factory	Model
Dehydrator	DIAPATH	Donatello
Embedding machine	Wuhan Junjie Electronics Co., Ltd.	JB-P5
Frozen platform	Wuhan Junjie Electronics Co., Ltd.	JB-L5
Constant temperature shaker	BLabotery	ZHPW-250
Electronic scale	Electronic	SF-400

1.2 Main experimental reagents

Reagent	Factory	Article Number
Xylene	Sinopharm GroupChemical Reagent Co., Ltd.	10023418
Absolute ethanol	Sinopharm GroupChemical Reagent Co., Ltd.	100092683
EDTA decalcification fluid	Servicebio	G1105

2. Experimental steps

2.1 Material: place the printed labels and tissues one by one in a perforated PE tube or embedding frame, then put them in a bucket, pour the new decalcification solution, seal them, and put them in a constant temperature shaker. In the process, decalcification is carried out by constant temperature shaker decalcification. The replacement period of the decalcification solution is 2-3d.

2.2 Decalcification: Put the decalcification keg into the constant temperature shaker for decalcification, and the temperature is generally controlled at 25 degrees Celsius. Between 30 degrees Celsius. The rate of the constant temperature shaker is controlled at 110-120RPM, and the degree of decalcification is observed every two days. If the needle can be moved, the tissue is cut with a Lycra blade according to the material requirements, which can accelerate the softening speed.

2.3 Dehydrating wax: take out the softened bones, rinse with running water for an afternoon, put it into the dehydrating basket, and dehydrate in a gradient gradient alcohol in the dehydrator. Details as follow: 75% alcohol for 2h ; 85% alcohol for 2h; 90% alcohol for 1.5h; 95%

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Add: 5th Floor, C22 Building, Biopark, No. 388 Gaoxin 2nd Road, East Lake High-tech Developing Zone, Wuhan, Hubei, China, 430079 Website: <http://www.servicebio.com> Tel: 027-5111-3188

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Wuhan Servicebio Technology Co., Ltd.

alcohol for 2h; anhydrous ethanol I for 2h; anhydrous ethanol II for 2h; alcohol benzene for 40min; Xylene I for 40min; Xylene II for 40min; 65°C melted paraffin I for 0.5h at 65°C; Melted paraffin II for 1h; 65°C Melted paraffin III for 2h45min.

2.4 Burying: The wax-soaked tissue is embedded in the embedding machine. Put the melted wax into the box first, remove the tissue from the box before the wax solidifies and put it into the box according to the accordingly. Cooling at -20°C freezing table, the wax block is removed from the embedding frame and trimmed after solidification.

2.5 Sectioning: Place the trimmed wax block in a paraffin section, 4μm. thick Slice float on the spreader 40°C warm water will spread the tissue flat, slide will remove the tissue, 60°C oven baking sheet. The water is roasted with dry wax and stored at room temperature.