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Add: 5th Floor, C22 Building, Biopark, No. 388 Gaoxin 2nd Road, East Lake High-tech Developing Zone,
Wuhan, Hubei, China, 430079 Website: <http://www.servicebio.com> Tel: 027-5111-3188

Ammonium Ferric Sulfate Dyeing Experimental Report

1. Experimental equipment and reagents

1.1 Experimental equipment

Name	Manufacturer	Model
Frozen section machine	Thermo	CRYOSTAR NX50
Slicers	Shanghai Leica Instrument Co., Ltd	LEICA 819
Adhesion microscope slide	Servicebio	

2. Main experimental reagents

Reagent name	Manufacturer	Article number
Ammonium ferric sulfate	Sinopharm Group Chemical Reagent Co., Ltd	10001818
Glacial acetic acid	Sinopharm Group Chemical Reagent Co., Ltd	10000218
Concentrated sulfuric acid		

2. Reagent preparation

2.1 0.5% ammonium ferric sulfate: 2.5 g ammonium ferric sulfate dissolved in 100mL distilled water.

2.2 Glacial acetic acid and concentrated sulfuric acid mixture: glacial acetic acid: concentrated sulfuric acid = 1:1 .

3. Experimental steps

3.1 Ammonium ferric sulfate staining: the frozen sections were naturally dried, immersed in 2.5% ammonium ferric sulfate solution, and treated in a 37 °C water bath for 3 days. The slides were taken out and washed with distilled water for 3 times. The water around the tissues was dried with filter paper, and the tissues were naturally dried. The mixture of glacial acetic acid and concentrated sulfuric acid was dropped on the tissues, covered slides , and the excess liquid was wiped off.

3.2 Observed under microscope and took images.

4. The results were as follows

4.1 It is start to appear red or blue after dropping the mixture of glacial acetic acid and sulfuric acid. After about 1-5 minutes, the tissue gradually appears green, which indicates the presence of cholesterol and cholesterol ester. After about half an hour, the green gradually fades to brown.

5. Precautions

5.1 The mixture of acetic acid and sulfuric acid should be stored in refrigerator.

5.2 Do not Touch by hand in dyeing process.

5.3 During the dyeing process, pay attention to the change of tissue color on the slides.

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Wuhan Servicebio Technology Co., Ltd.