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Add: 5th Floor, C22 Building, Biopark, No. 388 Gaoxin 2nd Road, East Lake High-tech Developing Zone,
Wuhan, Hubei, China, 430079 Website: <http://www.servicebio.com> Tel: 027-5111-3188

抗荧光淬灭封片剂	servicebio	G1401
Tunel 试剂盒	Servicebio	G1502

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1.2 Major reagents

Name	Producer	Code
Ethanol	SCRC	100092183
Biodewax and Clear Solution	SERVICEBIO	G1128
PBS solution	Servicebio	G0002
Permeabilize solution	Servicebio	G1204
Proteinase K	Servicebio	G1205

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Add: 5th Floor, C22 Building, Biopark, No. 388 Gaoxin 2nd Road, East Lake High-tech Developing Zone, Wuhan, Hubei, China, 430079 Website: <http://www.servicebio.com> Tel: 027-5111-3188

DAPI	Servicebio	G1012
anti-fade mounting medium	Servicebio	G1401
Tunel assay kit	Servicebio	G1502

2. Procedure

2.1 Deparaffinize and rehydrate: incubate sections in 2 changes of xylene, 15-20 min each. Dehydrate in 2 changes of pure ethanol for 10 min each, followed by dehydrate in gradient ethanol of 95%, 90%, 80%, and 70% ethanol, respectively, 5 min each (extend deparaffinize time slightly in winter). Wash in distilled water.

2.2 Antigen retrieval: eliminate obvious liquid, mark the objective tissue with liquid blocker pen. Add proteinase K working solution to cover objectives and incubate at 37°C for 25 min. then wash three times with PBS (pH 7.4) in a Rocker device, 5 min each. (Method for configuring working solution of proteinase K, stock solution: PBS=1:9)

2.3 Permeabilization: eliminate excess liquid, add permeabilize working solution to cover objective tissue, then incubate at room temperature for 20 min. wash three times with PBS (pH 7.4) in a Rocker device, 5 min each. (The membrane breaking fluid is 0.1% triton. Configuration method, triton stock solution: PBS=1:1000)

2.4 Equilibrium at room temperature: After the slices are slightly dried, buffer is added to the tissues in the circle, and the buffer is incubated at room temperature for 10 minutes

2.5 Tunel reaction: Take appropriate amount of TDT enzyme, dUTP and buffer in the tunel kit according to the number of slices and tissue size and mix at 1:5:50 ratio. Prepare this reaction solution according to demand before use. Add this mixture to objective tissue placed in a flat wet box, incubate at 37°C for 2 h. be sure to keep the wet box moist by adding water.

2.6 DAPI counterstain in nucleus: wash three times with PBS (pH 7.4) in a Rocker device, 5 min each. Then incubate with DAPI solution at room temperature for 10 min, kept in dark place. .

2.7 Mount: wash three times with PBS (pH 7.4) in a Rocker device, 5 min each. Throw away liquid slightly, then coverslip with anti-fade mounting medium.

2.8 Photomicrograph: The sections were observed under a fluorescence microscope and images were collected. (DAPI ultraviolet excitation wavelength 330-380nm, emission wavelength 420nm, blue light emission; CY3 excitation wavelength 510-561nm, emission wavelength 590 nm, red light emission)

3. Results

The nuclei stained by DAPI are blue under the excitation of ultraviolet light, the tune3 kit is labeled with CY3 fluorescein, and the positive apoptotic nuclei are red