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Add: 5th Floor, C22 Building, Biopark, No. 388 Gaoxin 2nd Road, East Lake High-tech Developing Zone, Wuhan, Hubei, China, 430079 Website: <http://www.servicebio.com> Tel: 027-5111-3188

Cell (FITC) Tunel Assay Protocol (Cell Climbing Slides)

1. Apparatus and reagents

1.1 Apparatus

Name	Producer	Model
Glass microscope slide	Servicebio	
Vortex	Servicebio	MX-F
Micro-centrifuge	Servicebio	D1008E
Rocker	Servicebio	TSY-B
Pipettor	Dragon	KE0003087/KA0056573
Liquid blocker pen	Servicebio	G6100
Ortho-Fluorescent Microscopy	Nikon	NIKON ECLIPSE C1
Imaging system	Nikon	NIKON DS-U3

1.2 Major reagents

Name	Producer	Model
PBS solution	Servicebio	G0002
Permeabilize solution	Servicebio	G1204
DAPI	Servicebio	G1012
anti-fade mounting medium	Servicebio	G1401
Tunel assay kit	Servicebio	G1501

2. Procedure

2.1 Permeabilization: dry the cell climbing slides slightly. Mark the objective area with liquid blocker pen, where add 50-100 μ l of permeabilize working solution. Incubate for 20 min at room temperature. Wash three times with PBS solution, 5 min each.

2.2 Equilibrium at room temperature: After the slices are slightly dried, buffer is added to the tissues in the circle, and the buffer is incubated at room temperature for 10 minutes

2.3 Tunel reaction: Take appropriate amount of TDT enzyme, dUTP and buffer in the tunel kit according to the number of slices and tissue size and mix at 1:5:50 ratio. Prepare this reaction

solution according to demand before use. Add this mixture to objective tissue placed in a flat wet box, incubate at 37°C for 2 h. be sure to keep the wet box moist by adding water

2.4 DAPI counterstain in nucleus: wash three times with PBS (pH 7.4) in a Rocker device, 5 min each. Then incubate with DAPI solution at room temperature for 10 min, kept in dark place.

2.5 Mount: wash three times with PBS (pH 7.4) in a Rocker device, 5 min each. Throw away liquid slightly, then coverslip with anti-fade mounting medium.

2.6 Microscopic examination and collecting images through fluorescence microscope. DAPI emits blue light at an ultraviolet excitation wavelength of 330-380 nm and an emission wavelength of 420 nm; FITC has an excitation wavelength of 465-495 nm and an emission wavelength of 515-555 nm, and emits green.

3. Results

Nucleus is blue by labeling with DAPI. Tunel assay kit is labelled with FITC. Positive apoptosis cells are green.

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Wuhan Servicebio Technology Co., Ltd.