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MMA Embedding

1. Experimental equipment and reagent

1.1 Equipment

Brand name	Manufactures	Model
Hard tissue slicer	Leica	HistoCore AUTOCUT
Disposable blades	Leica	TC-65
Water Bath	KEDEE	KD-P

1.2 Reagent

Reagent	Manufactures	Cat
4% PFA Fix Solution	Servicebio	G1101
70% FAA Fix Solution	Servicebio	G1103
Ethanol absolute	Sinopharm Chemical Reagent Co.,Ltd	100092683
Xylenes	Sinopharm Chemical Reagent Co.,Ltd	100023418
Methyl methacrylate	Macklin	M813511
Dibutyl phthalate	Sinopharm Chemical Reagent Co.,Ltd	10008518
Dibenzoyl peroxide	Macklin	B802244

2. Procedure for sectioning hard tissue

2.1. Draw materials

Fresh tissue was used to determine the sampling site and minimize mechanical damage such as pulling, contusing and squeezing. 4%PFA fixed bone tissue and 70%FAA fixed plants for at least 48 hours. It should be note that the tissue sampling thickness is about 5 mm and the maximum length is 3 cm. The bone tissue should try to shave the surface muscles, for fear that affect the observation of the results.

2.2. Dehydration

The tissue is then dehydrated in 60% - 70% - 80% - 90% - 95% alcohol. Each step takes anywhere for 24 to 72 hours. Next, absolute ethyl alcohol is used twice to dehydrate and 12 h to 24 h each time. The same is true for xylene.

2.3. MMA infiltration

Put tissue into three cylinders of permeates penetrating fluid at 4°C for 24 to 72 hours in turn.

2.4. MMA embedding

The tissue was put into the embedding bottle, and an appropriate amount of embedding agent was added. Then vacuumed for 5 h and polymerized in a 37°C water bath.

2.5. Sectioning

Routine sectioning about 10 μm by hard tissue slicer and place sections at 60°C oven overnight.