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1. Apparatus and Reagents

1.1 Apparatus

Equipment	Manufacturers	Model
Vertical refrigerated display cabinet	XingXing	LSC-316C
Horizontal freezer	MeiLing	BCD-318AT
Electronic balance	Mettler Toledo	ME203E/02
High speed tissue grinder	Servicebio	KZ-II
High speed refrigerated centrifuge	Dragon	D3024R
Electrothermal constant temperature water tank	Jinghong	DK-8B
Ultrapure Water Polishing System	Aiken water Elf	AK-RO-C2
Vortex mixer	Servicebio	MX-F
Enzyme label detector	BioTeK	Epoch
Automatic biochemical analyzer	Rayto	Chemray 240
Automatic biochemical analyzer	Rayto	Chemray 420
Automatic biochemical analyzer	Rayto	Chemray 800

1.2 Laboratory consumables

Equipment	Manufacturers	Model
1.5ml eppendorf tubes	Servicebio	EP-150-M
0-1000ul Adjustable-Volume Pipettor	Dragon	KE0037273
0-200ul Adjustable-Volume Pipettor	Dragon	YE3K030591
0-50ul Adjustable-Volume Pipettor	Dragon	DS35110
0-10ul Adjustable-Volume Pipettor	Dragon	KE0012951

2. Collection and preservation of samples

1. Serum samples (dry ice transport)

The blood sample should be placed at room temperature for 2 hours or 4 °C overnight, then turned at 2-8 °C 3000 RPM centrifuged for 15 minutes, and the supernatant can be detected immediately; or it should be repackaged and stored at - 20 °C or - 80 °C, but repeated freezing and thawing should be avoided. The thawed sample should be centrifuged again and tested.

2. Plasma sample (dry ice transport)

Heparin can be used as anticoagulant. Within 30 minutes after collection, the sample can be immediately detected at 2-8 °C for 3000 RPM centrifuged for 15 minutes; or it can be separately packed and stored at - 20 °C or - 80 °C, but repeated freezing and thawing should be avoided. The thawed sample should be centrifuged again and tested.

Note: the sample should be clear and transparent, and the suspension should be removed by centrifugation; hemolytic sample should be avoided. If the sample is not detected immediately after collection, it should be frozen in - 80 °C to avoid repeated freezing and thawing. There should be no bubbles in the sample during the test, so as to prevent the sampling probe from absorbing the bubbles and thinking that the sample has been inhaled. Hemolysis of the sample has an effect on the experiment. The sample size should be sufficient, and the detection needs to reserve a minimum amount (about 100ul).

3. Composition and storage conditions of kit

1. Main components: R1, R2

2. Storage conditions and validity period: the original kit is stored at 2-8 °C, valid for 12 months

4. Reagent preparation

Single reagent method is to mix R1 and R2 according to the proportion in the instruction manual or single reagent is to be used directly; double reagent method is to apply R1 and R2 respectively.

5. Operation steps

Kit: Rayto / Changchun Huili

5.1 Sample receiving: serum, plasma

5.2 Experimental steps:

5.2.1 Preparation of working reagent: prepare and mix R1 and R2 according to the proportion in the instructions by single reagent method; apply R1 and R2 respectively by double reagent method.

5.2.2 Set the corresponding parameters on the automatic biochemical instrument.

5.2.3 Sample Assay. Automatic biochemical analyzer.

5.3 Experimental results: the results are exported after the automatic biochemical analyzer is tested, and the results are sent in the form of Excel.

6. Normal reference range of rats and mice

Note: the reference value of serum biochemical index will be quite different due to the different conditions of animal breed, source, age, weight, feeding environment, nutritional status, inspection method, etc. The following reference values are calculated from the normal rat and mouse detection in the laboratory and are for reference only.

Rat	index	liver function					
		ALT	AST	ALB	ALP	GGT	TBA
	unit	U/L	U/L	g/L	U/L	U/L	umol/L
	Reference range	21.53-61.75	41.47-195.65	21.16-34.77	12.04-610.97	0.58-6.81	23.92-59.09
Rat	index	renal function			blood sugar	Other	
		BUN	CR	UA	GLU	TP	
	unit	mg/dL	umol/L	umol/L	mmol/L	g/L	
	Reference range	9.75-22.71	10.90-118.07	58.38-122.65	4.98-15.95	44.57-81.43	

Mouse	index	liver function					
		ALT	AST	D-BIL	T-BIL	ALB	ALP
	unit	U/L	U/L	umol/L	umol/L	g/L	U/L

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Add: 5th Floor, C22 Building, Biopark, No. 388 Gaoxin 2nd Road, East Lake High-tech Developing Zone, Wuhan, Hubei, China, 430079 Website: <http://www.servicebio.com> Tel: 027-5111-3188

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Wuhan Servicebio Technology Co., Ltd.